

THERMOCYCLER INSTRUCTIONS

1. Sign out thermocyclers PRIOR to using them.
2. Remove your name from the sign out book if you decide not to run a PCR.
3. Remove samples as soon as possible after your program has finished. Samples left in the thermocyclers will be placed in the refrigerator in Rm 231.
4. Do not use MJR thermocyclers for long incubations. Please use the Thermolyne thermocyclers for this.
5. Enter ALL programs onto a program log sheet.
6. Clean up any condensation that forms in cells when finished.
7. Keep thermocyclers in a staggered position to maximize air ventilation.
8. Manuals are located above the thermocyclers. Do not remove them from the lab.

Contact: Melanie Lewis Ivey (Rm 231, ext. 2704)